



RO System Flushing Instructions:

The manufacturer recommends draining RO system tank once per week through the RO faucet, and flushing 300ml of purified water through each dispenser daily. Failure to flush the system regularly in addition to intermittent or low volume system will compromise the quality of your system's dispensed product water.

If your system dispenses product water that contains mildew particulum, has a bad odor or taste, or contains any other type of suspended matter, then perform the following flushing procedure to renew the quality of purified water dispensed from your system.

To flush RO Tank & System:

- 1) Open RO faucet for 15 minutes to drain tank.
(RO faucet is the smaller drinking faucet installed on counter top).
- 2) Close RO faucet.
- 3) Wait 3-4 hours for tank to completely refill.
- 4) Dispense 600ml from each dispenser in pharmacy.
- 5) Open RO faucet for 30 minutes to drain remaining flushing water.
- 6) Wait 2 hours & evaluate new product water.
- 7) Repeat steps 1-6 if second flushing is necessary.
- 8) Flush minimum of 300ml through each dispenser daily.
- 9) Drain tank one time per week to ensure the highest purity of water.

~ RO system flushing procedure is now complete ~
If further assistance is required, call 1-800-464-8275.